

How to install and set up Repetier-Host (Bioprinting Assisting Software) on a Mac




Overview: The Repetier-Host Software allows Mac users to control their Bioprinters from their computers.

Materials:

You can find the link to download the latest Repetier-Host software [here](#)


Protocol:

1. Navigate to the Repetier-Host page by following the above link or locating the Repetier host link on the <https://cellink.com/support/> download page.

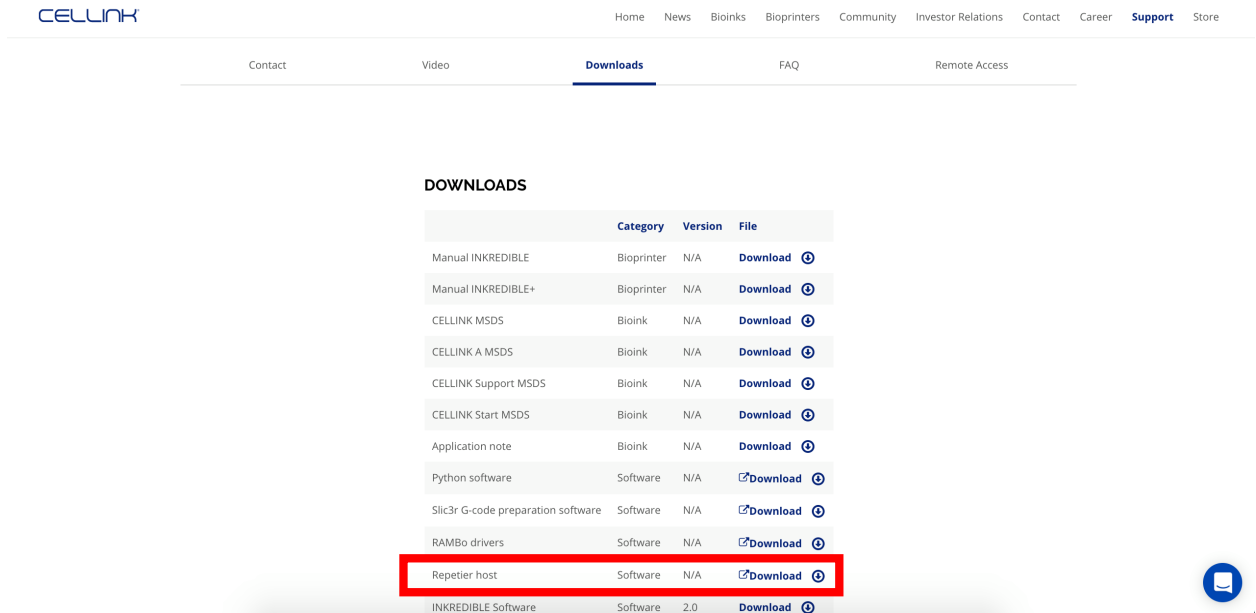

2. Download the version of Repetier-Host for Mac


3. Open the file and drag the icon to your Applications folder as shown in the photo

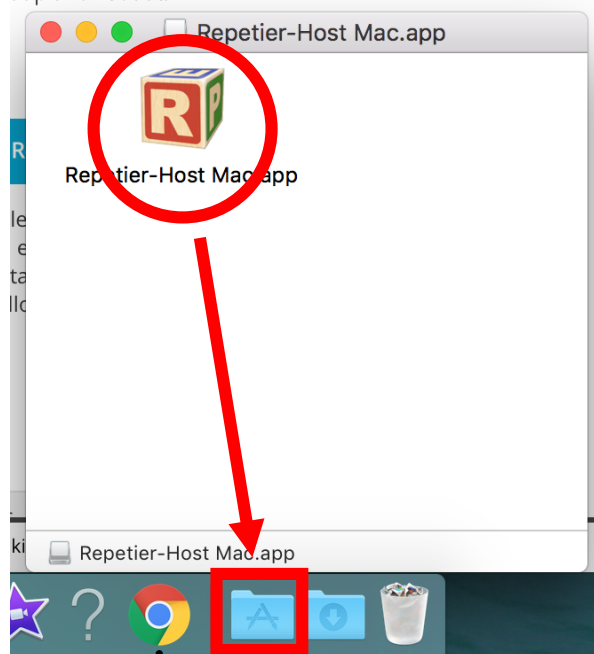


4. A security message may appear because the Repetier-Host Application is unfamiliar to the computer. To get around the security settings, open the application from the Applications folder in Finder

CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00

CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566

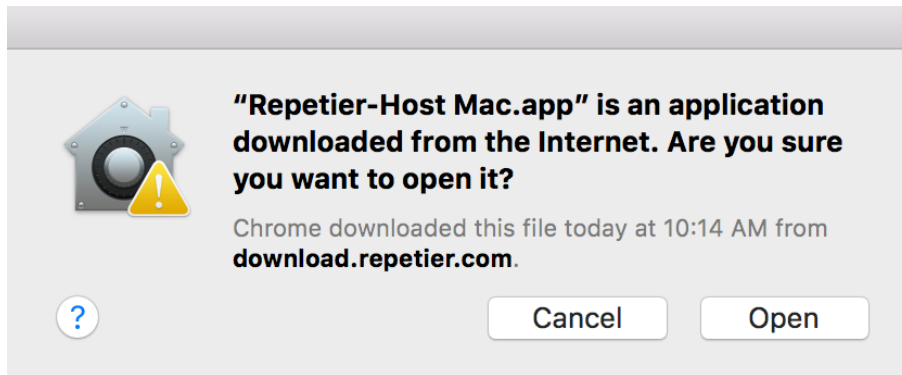


5. Click on "Printer Settings" and ensure that your printer is set to the following parameters

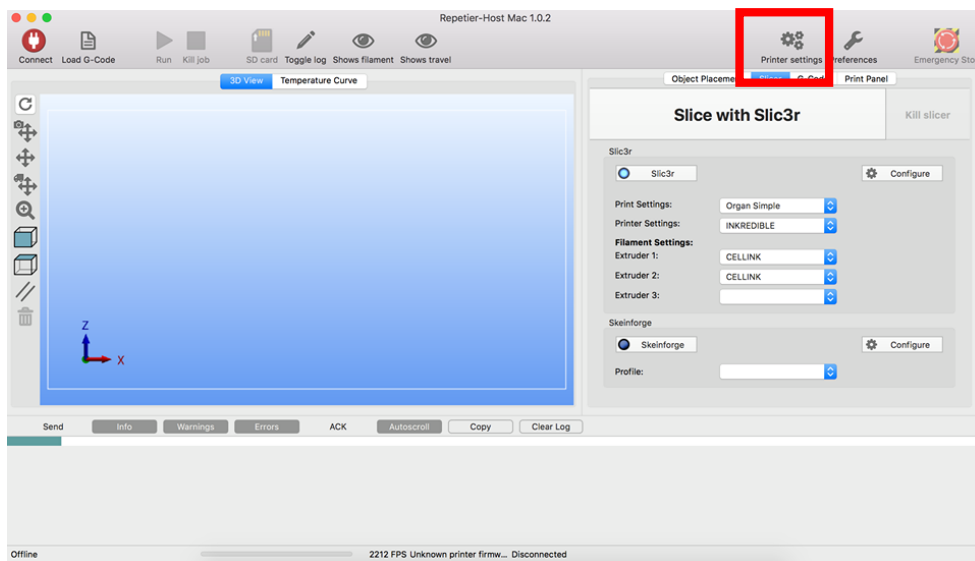



CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00

CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566

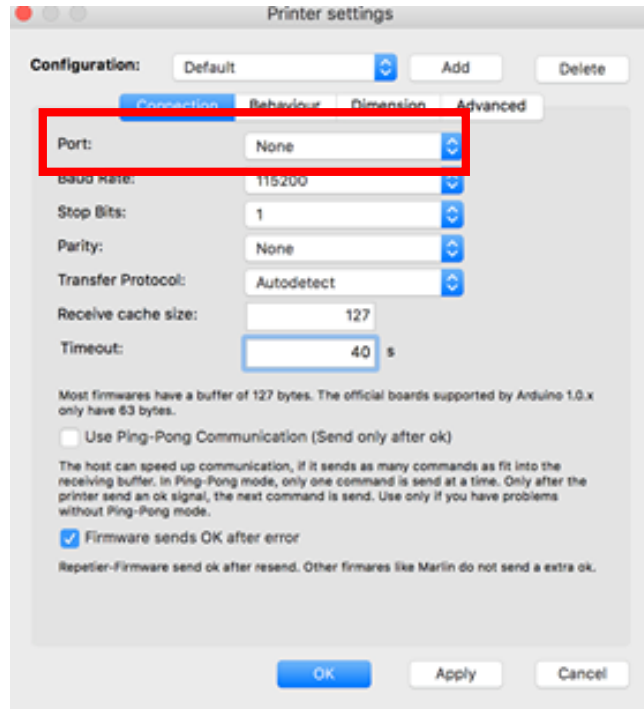


Under port, select the specific port that appears when the computer and printer are hooked up with the USB cable. For example “usbmodem1421”

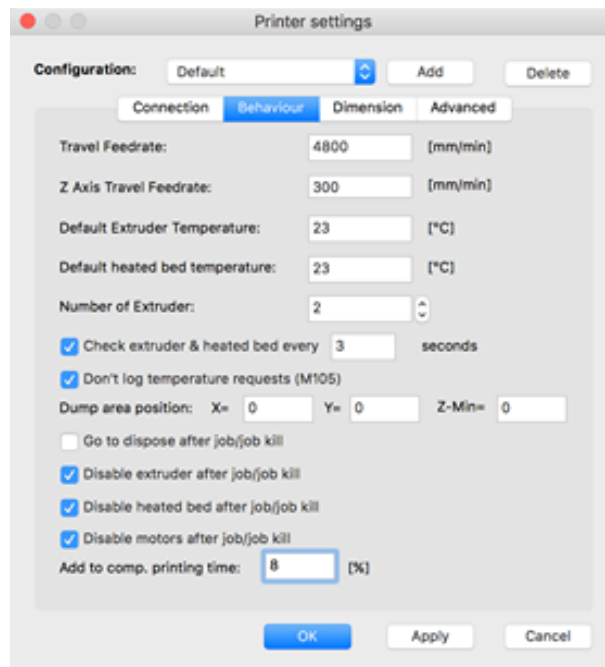

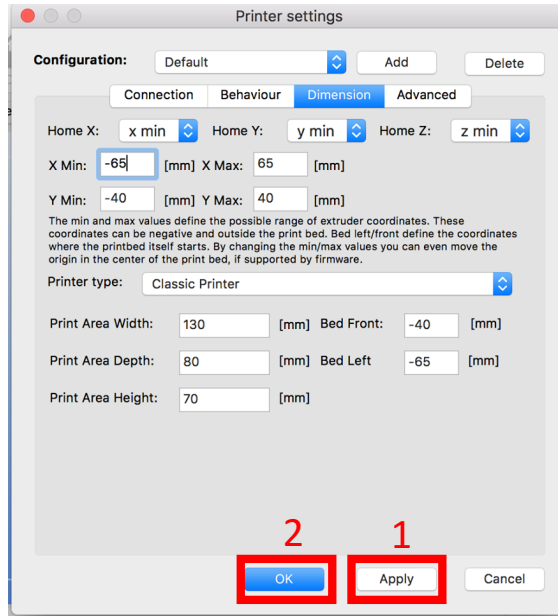

6. Once you are done, press Apply, then click OK
7. Ensure that the printer and the computer you are using are hooked up using a USB cable, then click the connect button on the top left.

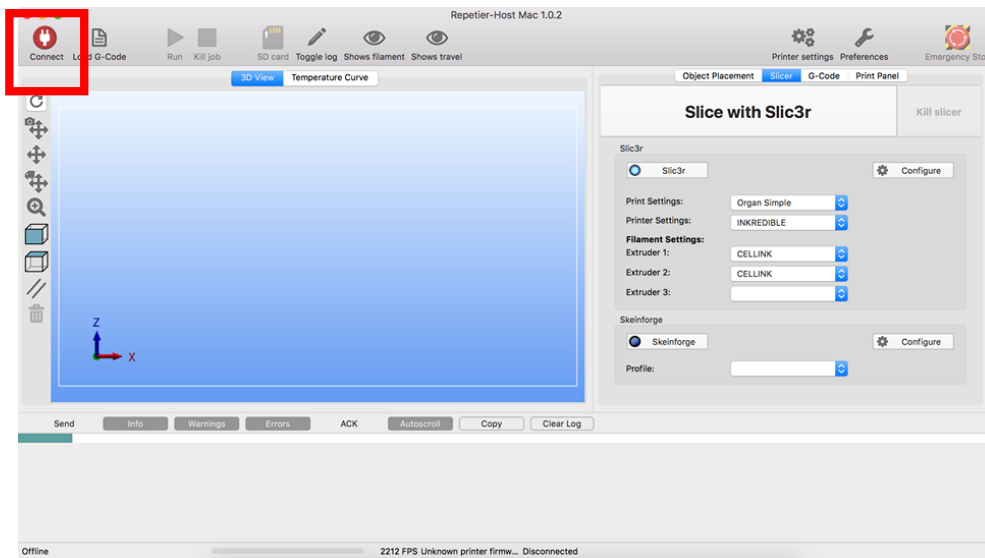

References:

Icon made by Freepik from www.flaticon.com

CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00

CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566

CELLINK®

Icon made by Smashicons from www.flaticon.com

Ref. No.:
Approved Date:
Version: 1

CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00

CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566